



Guia Rápido

Modelo de Vaca – PH42010101

O Modelo de Vaca é uma maquete realista em escala de um sexto, projetada para educadores da área agrícola ensinarem aos alunos sobre a anatomia bovina de forma prática e interativa. Este modelo realista apresenta características anatômicas internas e externas, incluindo os sistemas reprodutivo e digestivo, estruturas do úbere e muito mais.

Ideal para demonstrações em sala de aula ou estudo individual, o modelo permite que os alunos explorem a complexidade da anatomia da vaca de maneira envolvente e prática.



webASAP

R. Vinte e Oito de Setembro, 1413 – Vila
Dom Pedro I – São Paulo – SP, 04267-
000

Quer ter esse
**material
completo?**



[Fale com a gente
e saiba como!](#)



Guia Rápido

Modelo de Vaca - PH42010101



webASAP

R. Vinte e Oito de Setembro, 1413 - Vila
Dom Pedro I - São Paulo - SP, 04267-
000

Quer ter esse
**material
completo?**



Fale com a gente
e saiba como!



Guia Rápido

Modelo de Vaca – PH42010101



Small Cow Model Set

Quick Start Guide

The Small Cow Model Set includes five hands-on models for students to study the internal and external anatomy of a cow, including the ruminant, reproductive, and digestive systems.



Product includes:

- 5 Small Cow Models
- Full-color details
- 13 removable parts
- 1-year warranty
- Online access to curriculum, student activities and presentation slides



webASAP

R. Vinte e Oito de Setembro, 1413 – Vila
Dom Pedro I – São Paulo – SP, 04267-
000

Quer ter esse
**material
completo?**




**Fale com a gente
e saiba como!**

SMALL COW MODEL SET

Parts identification:

The items below make up the parts and pieces of each individual Small Cow Model.

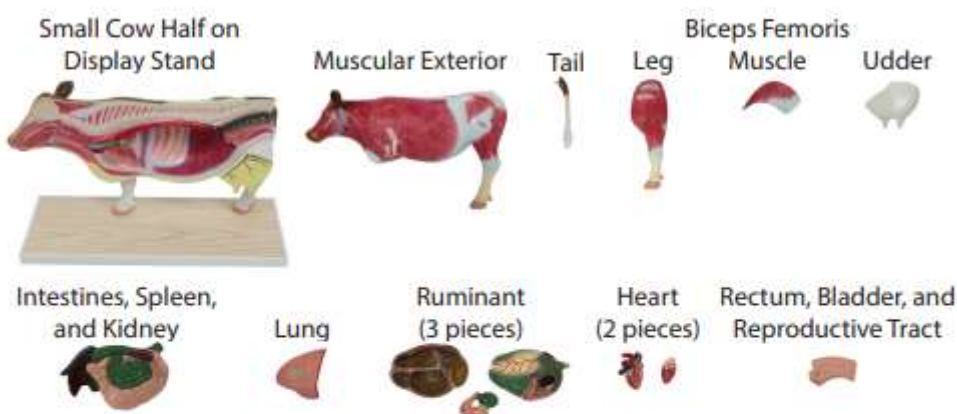

Assembly instructions:

Follow the instructions below for complete assembly of each Small Cow Model.

1. Begin with the small cow half on display stand
2. Insert the intestines, spleen, and kidney piece into the small cow half cavity
3. Connect the three ruminant pieces and insert them into the small cow half cavity on top of the intestines
4. Connect the two heart pieces and place them into the small cow half cavity
5. Place the lung onto the small cow half cavity, over the heart

